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Entlüftungsmodul HF.EA.8678

OPERATING INSTRUCTIONS

Extraction module HF.EA.8678

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Ontluchtingsmodule HF.EA.8678

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MANUAL DE INSTRUCCIONES

Módulo de ventilación HF.EA.8678

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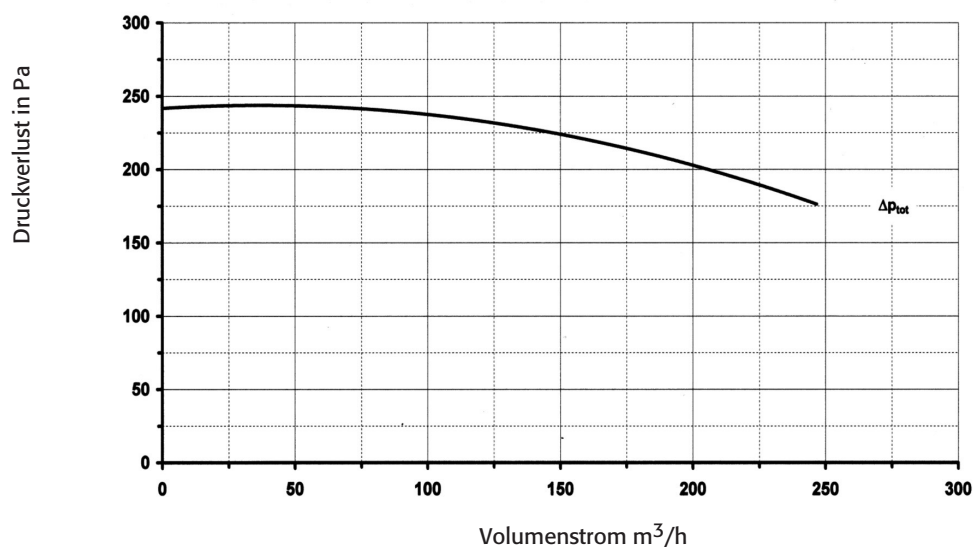
OPERATING INSTRUCTIONS

Dear Customer,

These operating instructions are for practical use and should be available to the user where the extraction module is to be used.

Keep these operating instructions best of all next to the extraction module.

An assured, faultless functioning of the extraction module is only ensured when these instructions are followed.

Do take note of the safety information.

Many thanks.

Your asecos Team

1. Development and manufacture

asecos GmbH Safety and Environmental Protection, D-63584 Gründau

2. Intended application

Extraction module for venting of up to four safety cabinets.

The device is used for the technical extraction of cabinets. The stream of exhaust air does not come into contact with electrically live parts or components with untreated metal surfaces. This means that the device is suitable for extracting fumes from cabinets which are defined in EN 14470 Parts 1 and 2. In addition, chemical cabinets as well as cabinets for storage of acids and alkalis can be vented.

3. Setting up the device

1. Bore two holes horizontally next to one another above the safety cabinet to be vented with a suitable drill at a distance of 500 mm (Ø 8 mm) from one another.
2. Blow any residues out of the drill holes and insert the two universal plugs supplied.
3. Screw the two screws supplied into the plastic plugs with a Phillips screwdriver until the screw heads are approx. 7 mm proud of the wall surface.
4. Hang the device onto the screw heads into the keyholes provided. Make sure that the shaft of each screw slides into the slot of the keyhole.

4. Connect device

1. Connect the suction holes (Ø 100 mm) to the 100 mm tubes on the suction nozzles of the safety cabinet. A reducer should be used for direct connection to the cabinet if necessary.
2. If you do not need one of the two 100 mm suction holes, please fit a blind plug (not included in the scope of delivery).
3. Connect the venting opening (Ø 140 mm) with the ventilation ducts on site.

5. Commissioning the device

1. Insert the mains plug into a socket that is earthed in accordance with regulations. (Note: The power supply cable attached to the device is 4 m long.)The green power supply control lamp goes on.
2. Operate the key switch in order to start the motor. The green operating lamp goes on. The monitoring electronics are activated at the same time.

6. Function control

The green power control diode must light up if everything has been correctly installed. The red monitoring control diode only lights up if the extraction volume is not correctly adjusted (e.g. because the extraction flap is closed or one or several cabinets are not exhibiting the necessary ventilation parameters). An acoustic signal is given at the same time. Acknowledge the signal and eliminate the cause .

7. Correcting faults if the extraction monitoring gives an alarm

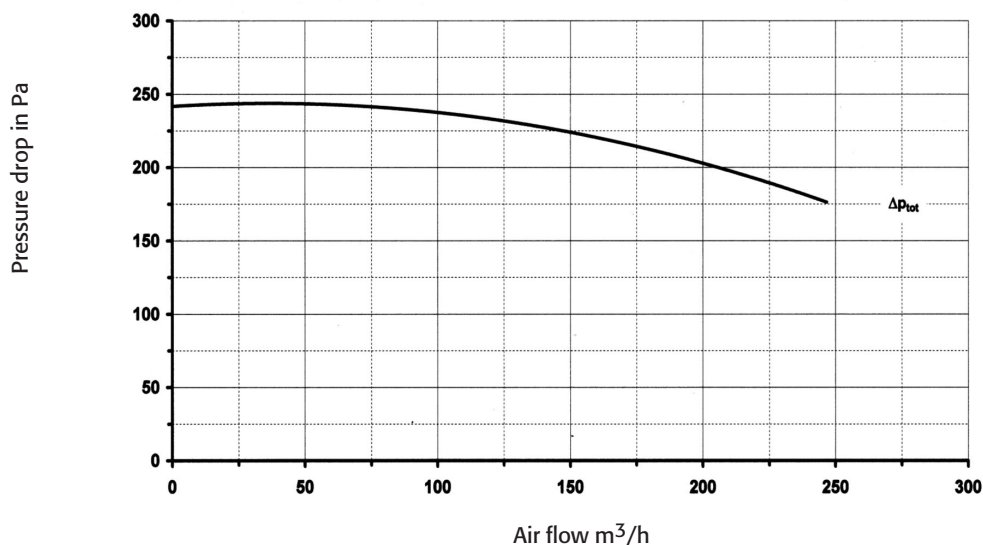
We can only take over a guarantee for perfect functioning of the device if it was correctly installed by a specialist firm authorised by us for the purpose.

Extraction monitoring works with a battery buffer, i.e. it will also give the alarm if the power supply is cut off. The alarm is given both visually and acoustically if the pre-set extraction parameters are changed by more than 10 %. After the alarm has sounded you can silence it initially by acknowledging the alarm (push reset). Then proceed as follows to locate the fault:

1. Check if the green LED (mains supply) is illuminated. If so, continue with 2.If not, check if the device is plugged into the mains socket, and if the socket is live.
2. Using a smoke bomb or other type of tracing vapour, check if air is being transported from each cabinet to the extraction nozzle of the device. The extraction hose provided on site has to be removed for this purpose. Fill the interior of each cabinet with vapour and look to see if the vapour becomes visible at the end of the device. If so, continue with 3.If not, the air passage is blocked. Please consult your ventilation service provider.
3. Have the volume of air which is extracted measured by the ventilation service provider. If the volume is correct, the pressure monitoring instrument (manometer) of the device should be adjusted. If the measured value does not correspond to the specifications, the air hose should be checked for foreign bodies and leaks. Remove all foreign bodies and seal all leaks which self-adhesive woven tape and close any remaining holes or joints with an elastic sealant. Start again with 1.If the fault is still present when you reach 3, please inform a specialist ventilation engineer or company.

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