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OPERATING INSTRUCTIONS

Dear customer,

These operating instructions are for practical use and should be available to the user where the suction hood is to be used. Please keep these operating instructions close to the unit. An assured, faultless functioning of the suction hood is only ensured when these instructions are followed.

Do take note of the safety information.

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2. Serial number

The serial number can be taken from the type plate on the front side of the suction hood.

3. Intended purpose

Complete capture of hazardous vapours, gases or suspended solids at the place where they emerge or where they are created before they can have effects that are damaging to health or to the environment (see hazardous materials regulations, workplace regulations and laboratory guidelines). Proof by means of ventilation test according to EN 14175-3 (5.4.4)

4. General safety notes / guarantee

- Take notice of the legislation and provisions to be applied in dealing with hazardous materials.
- Use the suction unit only if it is in good condition.
- Users must be instructed how to use the suction unit.
- Before using aggressive substances, please check the resistance of the surface of the suction unit to these substances.
- Any hazardous materials that escape must immediately be collected and removed.
- Take note of all information in these operating instructions.
- Please follow the instructions of the safety specialists, technical monitoring services (BG, GAA), construction supervision etc.
- Please observe accident prevention and workplace regulations.
- The required exhaust air flow is to be provided on site.
- **The guarantee for this product is agreed between you (the customer) and your dealer (the seller). Independently of this, the manufacturer provides a guarantee for the products listed in the operating instructions for 24 months after the date of sale. All model safety equipment are subject to a compulsory annual inspection by specialised staff authorised by the manufacturer. If this inspection is not carried out annually, or not carried out by specialised staff authorised by the manufacturer, the customer's guarantee claims against the manufacturer will be invalidated.**

5. Design

- Anodised aluminium tubular frame design with painted aluminium corner connectors
- The sides, ceiling, front and rear wall are made of transparent 5 mm-thick acrylic glass.

6. Function

Due to extraction at the ceiling, the suction unit is a highly effective way of ensuring that no vapours, gases or suspended solids involved in handling hazardous work materials (for instance when filling containers, gluing, cleaning, preparing, weighing etc.) pollute the breathable air. These are collected by the effective suction and removed by the exhaust system.

7. Commissioning

Testing before setup:

An appropriate exhaust system must be connected on site to the outlet connector (DN160) for the extraction of hazardous substances. Please refer to section 10 "Technical data" for the ventilation values

Commissioning:

- The ventilation values (see section 10 "Technical data") of the suction unit have been determined by the ventilation test method according to EN 14175-3 (5.4.4).
- Establish an on-site extraction air connection. A negative influence on the flow behaviour in the suction unit must be avoided. Depending on the place of installation and the possible disturbances associated with it, considerably larger extraction air volumes may be necessary in practice than the values set during the ventilation test.
- Open windows and doors at the place of installation (creation of air currents)
- Incorrect operation (i.e. no fast movements > 1 m/s)

8. Maintenance, care

Cleaning with normal mild household cleaners. Escaped fluids should be collected / removed immediately.

9. Ventilation adjustments

Efficient, secure retention of the pollutants in the suction unit is achieved through an optimum combination of fresh and extraction air.

Please note:

- The retention capacity of the suction unit depends on the air speeds achieved for the fresh air stream of the device.
- The determination of optimised fresh and extraction air flow ratios is significantly influenced by air movements present at the place of installation.

10. Technical data

The figures in the table are provided as a guide, and are not to be used in the sense of assured properties.

Model	ASE.50.6.4	ASE.50.9.4	ASE.60.9.6	ASE.90.12.6
External dimensions in mm:				
Height	500	500	900	900
Width	600	900	900	1200
Depth	400	400	600	600
Front opening in mm:				
Width	484	600	600	900
Height	437	437	542	542
Work surface in mm:				
Width	567	867	867	1167
Depth	383	383	583	583
Weight (kg)	11,6	14,6	26,4	30,4
Extraction air connection (NW/mm)	160	160	160	160
Recommended extraction air flow (m ³ /h)	250	300	360	540
Pressure loss (approx Pa)	45	64	93	208

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